

#### Shutdown protocol (Tubes)

1. Set sample flow switch to Hi
2. Add 2mL of 10% bleach to a test tube (do not fill more than  $\frac{1}{2}$  way)
3. Load tube and run 5 minutes
4. Add 2mL DI H<sub>2</sub>O to a test tube (do not fill more than  $\frac{1}{2}$  way)
5. Load tube and run for 5 minutes
- 6. Set fluidic switch to STNBY (standby)**
7. Leave the tube of DI H<sub>2</sub>O on the cytometer (except Canto)
8. If it after 5pm M-F, turn off

Turn off after all experiments Sat, Sun and on Holidays