

Exporting Bruker Data to ASCII files (20080108)

This guide covers exporting Bruker data from Topspin to ASCII files that can be plotted with programs like Excel and Origin.

To convert Bruker data to ASCII format:

1. Open Topspin and load the spectrum (file name, experiment #) you want to save.
2. Type 'convbin2asc'. This will start a program that will convert your spectrum. Once it is finished, close TopSpin.
3. Go to My Computer and then to Z:/data. Open the folder for your <userID>/nmr/<file name>/<experiment #>/pdata/1. Within that last folder, should be an ASCII file named 'ascii-spec'.

You could also open your Bruker spectrum in MestRec and just export it as ASCII. Might be simpler that way.