

## Exporting Bruker Data to Graphic files (20080108)

**This guide covers exporting Bruker data from Topspin and Bruker's layout software, Topspin Plot Editor on the PC computers in the off-line processing room.**

Both Topspin and the Topspin Plot Editor can create graphic image files from the NMR data displayed in the program.

The DCIF recommends using the saving your graphic files on the network drive Z:  
Z:\data\UID\filename.ext

Scalable Graphics such as .ps, .pdf, .wmf, .emf do not loose resolution when resized.

Pixilated Graphics such as jpg, .pcx, .png, .tif, .bmp may loose resolution when resized.  
.png are recommended for web publishing ( the use of jpg is not recommended)

### **TOPSPIN**

1. Arrange the spectrum as desired
2. Select File > Export
3. Type in the filename **and** the desired extension (file type)  
Example:  
filename.tif
4. Click [OK]

The possible extensions are: .png, .jpg, .jpeg, .bmp, .emf, .wmf

### **TOPSPIN PLOT EDITOR**

1. Arrange the spectrum as desired
2. Select File > Export
3. Type in the filename and select the desired extension (file type)  
The program will append the desired extension to the filename
4. Click [OK]

The possible extensions are:.emf, .ps, .pdf, .jpg, .pcx, .png, .tif