

Opening Varian data in MestReC

1. Using SecureFX, transfer in binary mode the data folder (ends in *.fid*) from Zippy to your folder (user ID) on the Z:\ drive, or to whatever other destination you prefer.
2. Remember that each Varian *name.fid* folder contains 4 subfiles: *fid*, *procpa*, *log*, and *text*.
3. Open MestReC.
4. Go to **File** → **Import Spectra**.
5. Find and open the *name.fid* folder.
6. Select the *fid* file within that folder, and click Open.

Opening Bruker data in MestReC

1. Using SecureFX, transfer in binary mode the data folder from the B400, B401, or B600 computer to your folder (user ID) on the Z:\ drive, or to whatever other destination you prefer.
2. Open MestReC.
3. Go to **File** → **Import Spectra**
4. Find your data folder.
5. When you originally analyzed your sample, you had assigned a name and an experiment number to each spectrum you collected. Within the data folder, you will now see the experiment number folders.
6. Open the experiment number folder that contains the spectrum you want to process.
7. Select the *fid* file and click Open.